Fluid bed systems



We set the standard

The very best for your laboratory

The innovative GPCG 2 LabSystem from Glatt

Absolute flexibility

The full range of fluid bed processes. In top quality. For new ideas and top products.

- Drying (in the granulating insert).
- Granulating (Top Spray in the granulating insert).
- Coating (Bottom Spray in the HS-Wurster*).
- Pelletizing (Tangential Spray in the rotor insert).

Optimum batch size - from 0.3 kg to 5 kg**. With textile filter or metal filter and various bottom screens. With Glatt single-arm spray nozzle, powder-feed and other innovations. In the standard version or customised to suit your specific requirements. The perfect equipment to ensure your success.

Step up safety

Your safety is of utmost importance to us. The PRO concept from Glatt. No compromises instead maximum safety.

- No pressure relief.
- Environmental and personal protection with a contained system.
- For all dusts up to an explosion pressure of 10 bar and their hybrid mixtures.
- · New sealing system. Less weight and better handling.
- As a total containment version for highly active substances.

Top for clinical sample production

The highest of requirements apply in the manufacture of clinical samples. Problems are something you can do without. What you need is the GPCG 2.

- Easy-to-clean design.
- Comprehensive process documentation.
- · Validation of processes.



Batch sizes from 0.3 kg to 5 kg**









Function and form taken to perfection

The innovative GPCG 2 LabSystem from Glatt

Optimum process control

Customer expectations are high. Fast and simple process control is what counts. The solution: ProcessView from Glatt.

- Touch-screen control.
- · Swivelling operating panel.
- Perfect ergonomics for the operator.
- Integrated recipe handling.
- Batch report.

Excellent handling and integrated technology

Perfect solutions and sophisticated design. The GPCG 2 offers all the advantages of integrated technology and top quality processing.

- Centralised process inserts, removeable and rotatable for product discharging.
- Simple filling and discharging.
- · Problem-free cleaning.
- Plug & Play. Just connect it and you're off.
- · Ergonomically arranged connections.
- Rotating mesh for contained emptying. Removable product container.
- Unique work surface made of synthetic granite. Resistant to solvents, scratch-proof and impact-proof. It still looks like new even after years of use.

Modular design

Expect absolute flexibility. Integrated modules offer you options now or in the future.

- Dehumidification of inlet air.
- Downstream deduster for dust sensitive coating processes.
- · Additional storage space for process accessories.





Swivelling operating panel







Removeable and rotatable product container


Technical Data / Addresses

GPCG 2 LabSystem

Batch sizes		small 1)	standard	large	
	Granulating insert	1 - 3 I	2 - 61	5 - 12	
	Wurster insert (4" and 6") 2)	21	3.5	_	
	Rotor insert 2)	-	51	-	
Electrical power ³⁾	Operating voltage	1 x 230 V and 3 x 400 V			
	Frequency	50 Hz			
	Total connected load		3.5 / 15 kW		
	Connector	CEE	CEE form 3-pin / 5-pin		
Compressed air	Max. consumption		12 m³/ h		
	Pressure		6 bar(g)		
	Connection	Н	Hose connection		
Inlet air heater, electric	Max. inlet air temperature	120° C			
Dimensions	Height		2427 mm		
	Width		1900 mm		
	Depth	987 mm			
Weight		1000 kg			
Spray pump 2/3	Max. output		53 ml / min		
	Connection voltage		230 V		
	Frequency		50 Hz		

¹⁾ Only together with the tower bypass for process air 2) Optional 3) Other voltages and frequencies on request

(Subject to technical alternations)

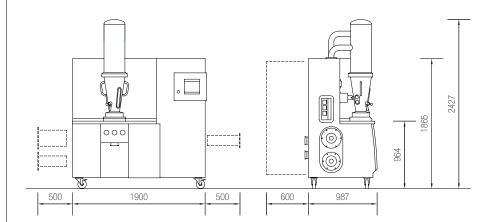

The innovative GPCG 2 LabSystem from Glatt

- Drying
- Granulating
- Coating
- Pelletizing





Glatt GmbH Process Technology

Werner-Glatt-Straße 1 79589 Binzen / Germany Phone: +49 7621 6 64 0 Fax: +49 7621 6 47 23 eMail: info@glatt.com

Glatt Maschinen- & Apparatebau AG

Kraftwerkstrasse 6 4133 Pratteln / Switzerland Phone: +41 61 8 26 47 47 Fax: +41 61 8 26 48 48 eMail: info@glatt.com

Glatt Ingenieurtechnik GmbH

Nordstraße 12 99427 Weimar / Germany Phone: +49 3643 47 0 Fax: +49 3643 47 12 31 eMail: info@glatt-weimar.de

Glatt Air Techniques Inc.

20 Spear Road Ramsey, NJ 07446 / USA Phone: +1 201 8 25 87 00 Fax: +1 201 8 25 03 89 eMail: info@glattair.com

Glatt Norden ApS.

Skøjtevej 27 - 31 2770 Kastrup / Denmark Phone: +45 48 14 22 44 Fax: +45 48 14 22 55 eMail: info@glattnorden.dk

Glatt Pharmatech SASU

Parc Technologique 6, rue Louis Néel 21000 Dijon / France Phone: +33 3 80 74 32 64 Fax: +33 3 80 74 43 95 eMail: info@glatt-pharmatech.fr

Glatt Protech Ltd.

Swannington Road
Cottage Lane Ind. Est.
Broughton Astley
Leicester LE9 6TU / Great Britain
Phone: +44 1455 28 58 58
Fax: +44 1455 28 55 10
eMail: info@glatt-protech.co.uk

www.glatt.com

